


Note: It may be necessary to press down on the cover through the slot.



2. Uncap the reagent bottle and screw it into the dispenser.

The reagent is now considered open for calculating the expiration time.

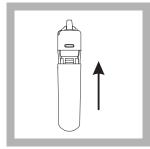

3. Replace the dispenser cover by sliding it into the grooves on the body until it snaps into place.




4. To dispense the powdered reagent, invert and grip so that the index finger is over the trigger. Engage the trigger fully, then release it quickly.


The blue indicator button is flush with the dispenser body when the trigger is completely released.



IN THE U.S.A., CALL TOLL-FREE 800-227-4224 FOR TECHNICAL INFORMATION AND ORDERING

Maintaining the SwifTest™ Dispenser

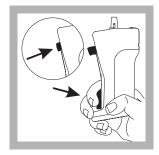

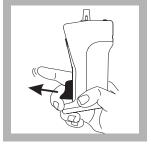





1. If the blue indicator button doesn't retract when the trigger is released, hold the dispenser with the dispensing tip up and engage the trigger fully.

2. Slide the finger off of the trigger so that it releases with a "snap". Snap the trigger until the indicator button retracts fully when the trigger is released. Invert to continue dispensing the reagent.

Note: Releasing the trigger quickly during reagent dispensing helps prevent powder build-up.

3. When the reagent bottle is empty, remove the dispenser cover and unscrew the empty bottle. Shake the excess powder out, then replace the dispenser cover. Snap the trigger 2 or 3 times over a disposal bin or lab sink.

Note: Read the reagent MSDS for disposal instructions.